

ez-Migration Trap (DC0019)

MegaView Science Co., Ltd.


<http://MegaView.com.tw> • <http://BugDorm.com>

Package Contents

- 1 ez-Migration Trap
- 2 Support Poles
- 2 Bottle Connecting Rings
- 6 Collecting Bottles
- 8 Guy Ropes
- 14 Tent Pegs
- 1 Carrying Bag

Assembly

1. Assemble poles. Make sure ends are completely seated in joints.
2. Push ends of one pole into pockets of one end of ez-Migration trap.
Careful: do not over bend pole.
3. Clip hooks around pole. Take care to maintain gentle curve in pole.
4. Repeat Steps 2-3 with the other pole and the other end of ez-Migration trap.
5. Stake out guy ropes from center of each end (where black changes to white) of ez-Migration trap to make it stand.
6. Stake base of ez-Migration trap, especially at ends of poles.
7. Remove moth excluder from collecting head entrance if not needed.
8. Additional guy points are sewn into ez-Migration trap for use in wind.

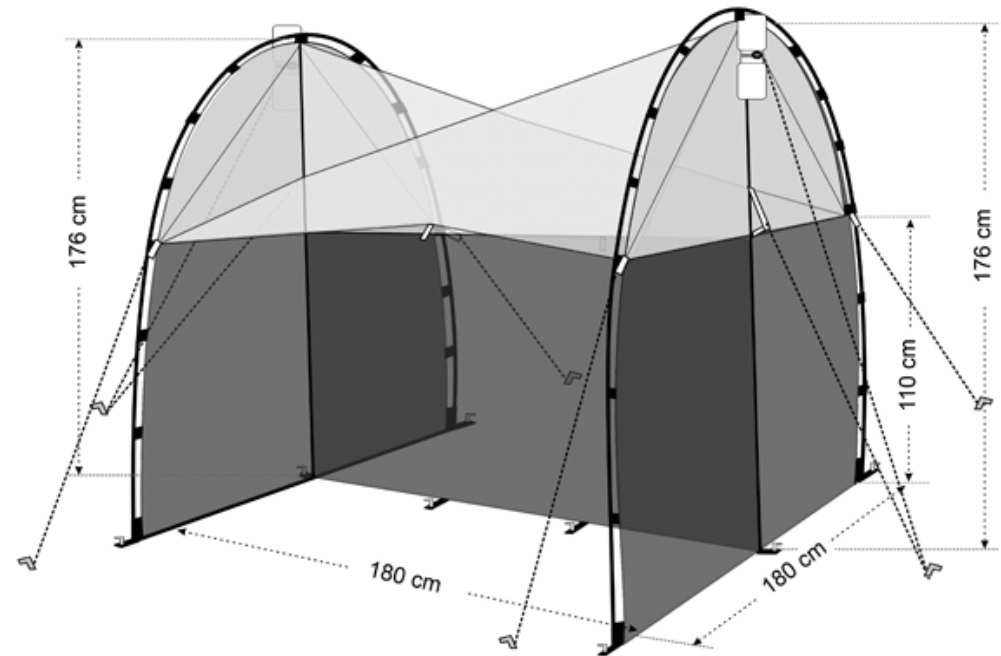

Have any questions? Please email us: bugdorms@megaview.com.tw